

DRIVER & SOFTWARE INSTALLATION

■ Driver


1. Connect the camera with PC via USB port, “Found New Hardware Wizard” appears. Choose the first option **“Yes, this time only”**, then click **“Next”**



2. Choose **“Install from a list or specific location (Advanced)”**, click **“Next”**.



3. Insert Driver/Software CD, check **“Include this location in the search”**. Browse the content on CD and select the directory **“H:\Driver_Plug_Play”** (where H denotes the drive letter in use). Click **“Next”** to continue.

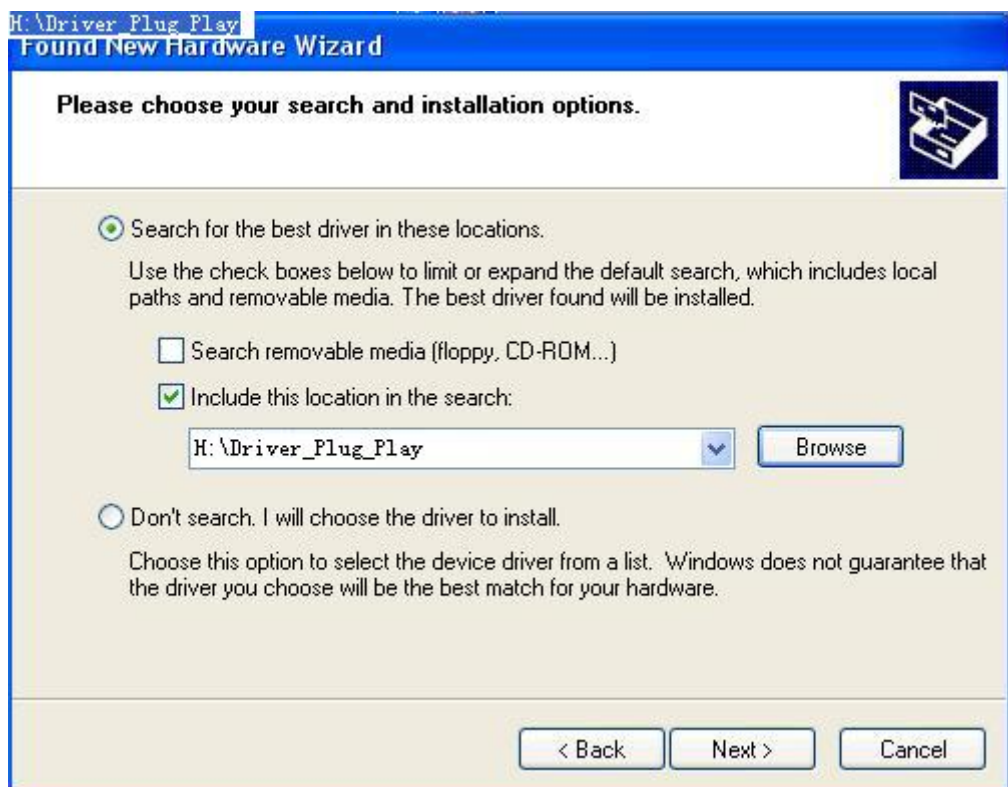

4. Click “Continue Anyway” when the following message prompts.



5. Click “Finish” to complete the installation process. New hardware device “USB2.0-CAM” has been successfully installed.




■ Software

1. Double click “**setup.exe**” to begin installation. Click “**Next**”.



2. Specify a folder for installing the software. Default path is “**C:\Program Files\ScopeView**”. Then click “**Install**”.

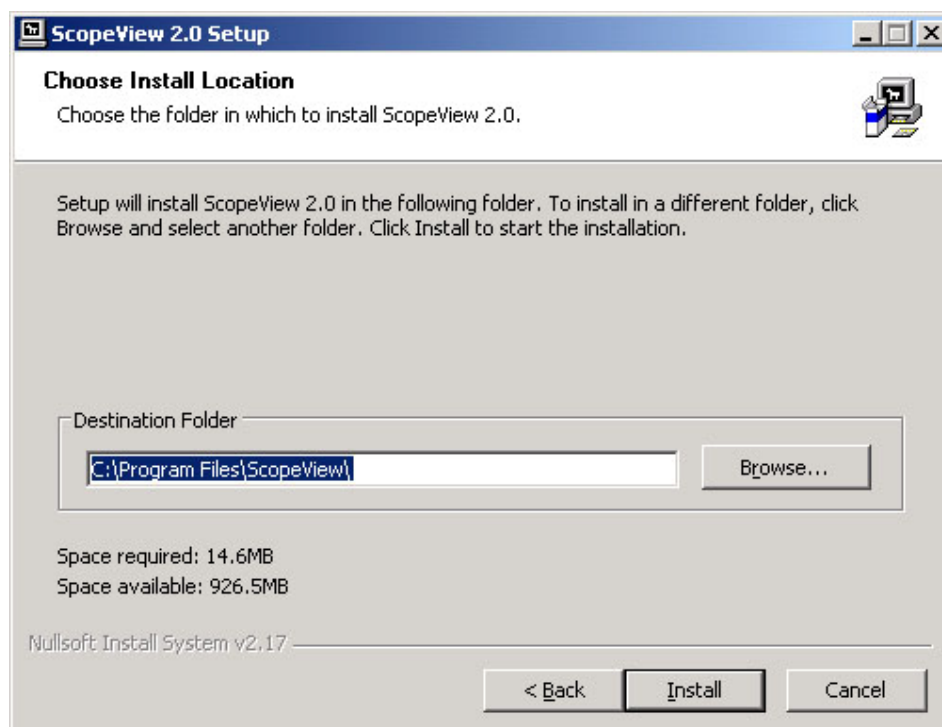

3. Click **“Finish”**. Installation is now done. Check **“Run ScopeView 2.0”** to launch the program upon exit.

